



## Event Calendar Add-on



API manual

2021/07/08

# Summary

Create calendar.....	3
Get calendar.....	3
Create event.....	3
Add external links to event .....	4

# Create calendar

**POST** http://[host]/extra\_modules/eventcal/API2/calendar.api.php

Headers	
X-LC-APP-Auth *	[The token generated in LC API setup]

  

Data	
actiontype *	createCalendar
uid *	[Contact ID] ex.: "1"
name *	[Calendar name] ex.: "My Calendar"
access_contacts	[Contact IDS separated with comma] ex.: "2,6,7", or "2"

**Returns** calendar ID  
JSON "2"

# Get calendar

**GET** http://[host]/extra\_modules/eventcal/API2/calendar.api.php?id=[Calendar ID]

**Returns** calendar data  
JSON {"id":"1","user\_id":"1","name":"TestCal"}

# Create event

**POST** http://[host]/extra\_modules/eventcal/API2/calendar.api.php

Headers	
X-LC-APP-Auth *	[The token generated in LC API setup]

  

Data	
actiontype *	createEvent
cid *	[Contact ID] ex.: "1"
title *	[Calendar name] ex.: "My event"

start *	[Start date] ex.: "2021-08-01 09:30:00"
end *	[End date] ex.: "2021-08-01 09:30:00"
description	[Event description] ex.: "Description of my event"
recurring	[DAILY WEEKLY MONTHLY]
until *	[Until this date] ex.: "2021-08-15" (mandatory if until is specified)

**Returns** event ID  
JSON "18"

## Add external links to event

**PUT** http://[host]/extra\_modules/eventcal/API2/calendar.api.php

Headers	
X-LC-APP-Auth *	[The token generated in LC API setup]

Data	
actiontype *	addEventExternalLinks
event_id *	[Event ID] ex.: "1"
external_links *	[External links separated with comma] ex.: "extra_modules/stability_test_manager/index.php?action=view_plan&plan_id=16, extra_modules/stability_test_manager/index.php?action=view_plan&plan_id=17"