LabCollector

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End date

All day event No

Description Auto-generated with Stability Test Manager addon

Recurring Once

Color #0070C0

Linked URLs https://labcollector.online/agilebiol/extra_modules/stability_test_manager/index.php?action=view_plan&plan_id=6#action-view-test-112

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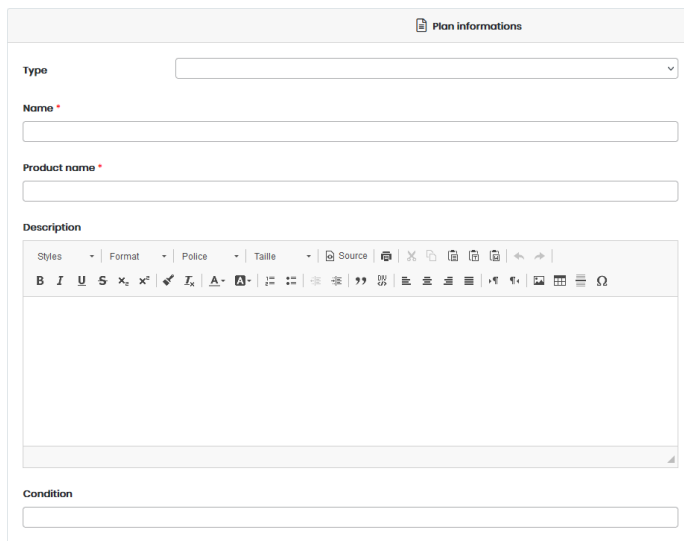
The screenshot shows a web form titled "Plan informations". It contains the following fields and elements:

- Type:** A dropdown menu.
- Nome ***: A text input field with a red asterisk indicating it is required.
- Product name ***: A text input field with a red asterisk indicating it is required.
- Description:** A rich text editor with a toolbar containing options for Styles, Format, Police, Taille, Source, and various text and image formatting tools.
- Condition:** A text input field.

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Start date * 2021-12-01 End date * 2022-12-31

1 Cycle * 1 Per * day AT * ex: 09:30
ex: 1 time a day at 9:30

2 Every * 30 days Per * 0 days
ex: every 30 days of 3 days

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1 Cycle * 1 Per * week AT * 15:30
ex: 1 time a day at 9:30

Choose 1 day(s) *

Monday Tuesday Wednesday Thursday Friday Saturday Sunday

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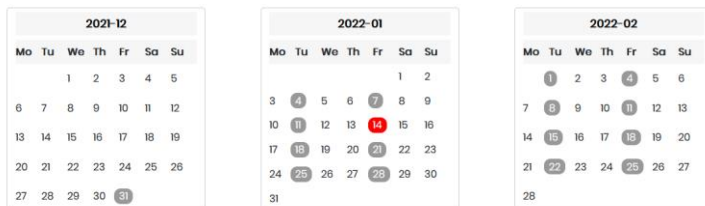
Planning preview

You can click on a date to exclude it from plan:

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The screenshot shows the 'Addons configuration' dialog for the Event Calendar module. It contains the following elements:

- Addons configuration** (header)
- Create a new calendar in Event Calendar addon** with radio buttons for Yes and No.
- Events color** with a text input field containing '#3788D8' and a color selection button (blue square).
- Connect to LSM** with radio buttons for Yes and No.

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
The screenshot shows the 'Addons configuration' dialog for the LSM module. It contains the following elements:

- Addons configuration** (header)
- Connect to LSM** with radio buttons for Yes and No.
- LSM URL domain *** with a text input field containing 'https://labcollector.online/agilebio1' and a sub-note: 'The first part of your LSM url, before extra_modules'. A green status indicator below reads 'LSM remote URL found'.
- LSM API token *** with a text input field containing '8790cd9b9aa4b888dea2d6aa2fed9db34918381c' and a sub-note: 'A token to allow LSM connection'.
- Group LSM orders by plan** with radio buttons for Yes and No.
- LSM tests** with a text input field and a sub-note: 'Enter at least 3 characters'. Below are three examples: '#2 Appearance x', '#3 Color x', and '#4 pH test x'.
- LSM sample type** with a text input field and a sub-note: 'Enter at least 3 characters'. Below is an example: '#2 Drug Powder x'.
- LSM model** with a text input field and a sub-note: 'Enter at least 3 characters'.

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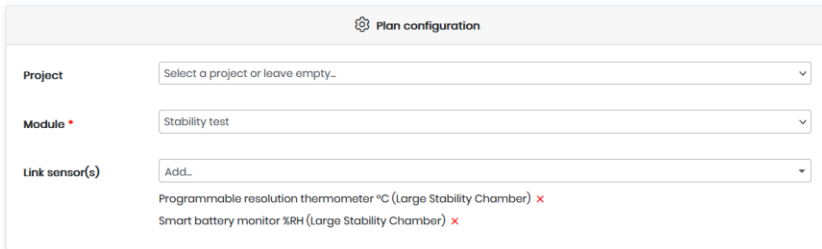4-2-3. Data logger

Plan configuration

Project: Select a project or leave empty...

Module: Stability test

Link sensor(s): Add...

Programmable resolution thermometer °C (Large Stability Chamber) x

Smart battery monitor %RH (Large Stability Chamber) x

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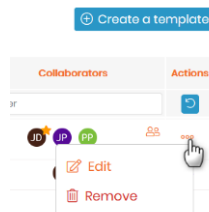
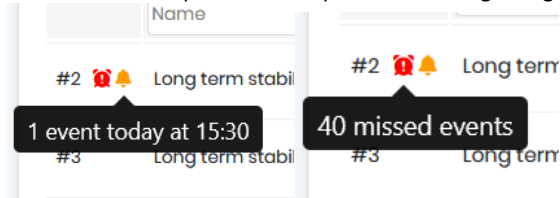
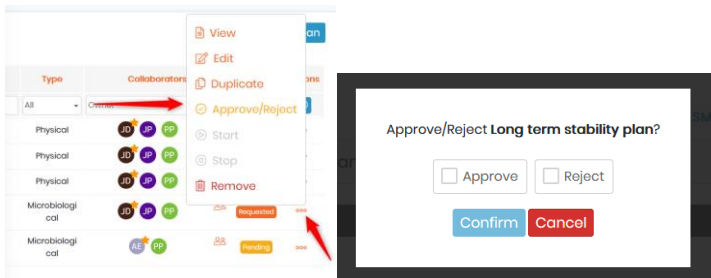
Box model > Storage location > Boxes prefix > Backup storage > Start

- 1 | Box 100 tubes (100 positions)
- 2 | Large stability chamber (+25°C), Rack1_6
- 3 | Long_term_plan_2022-1_
- 4 | Large stability chamber (+25°C), undefined

5 | **Confirm plan start:**

4 samples will be created in **Stability test** module.

2 backup samples will be created in **Stability test** module.

4 events will be created in **Long term stability plan** calendar (Event Calendar addon).

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Available events				
Missed events				
	Name	Product name	Available event	Action
#2	Long term stability plan	Product A	20/05/2022 event missed (late for 5 days)	
#2	Long term stability plan	Product A	17/05/2022 event missed (late for 8 days)	
#2	Long term stability plan	Product A	15/05/2022 event missed (late for 12 days)	


Do the sampling/test
Enter in the plan

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Long term stability plan: You have an event today at 15:30 (late for 2 hours and 12 minutes)

no	Start date	End date	Frequency	Project	Type

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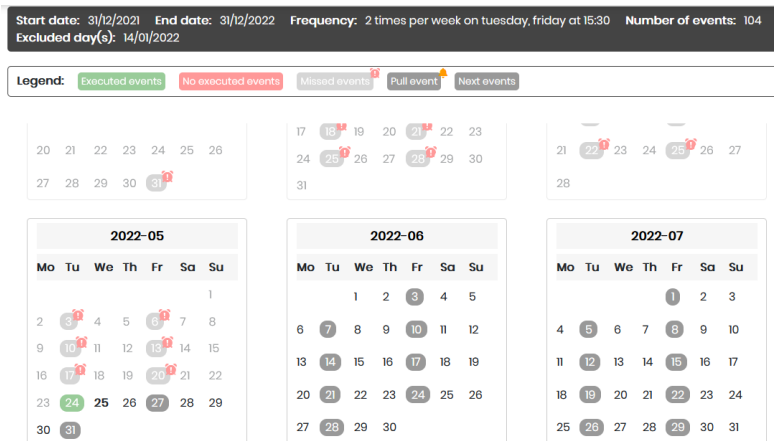113 SA_113

Stability Tests

Plan:	Short term stability plan
Test date:	23/05/2022
Status:	Executed (on 24/05/2022 18:02:54 by Jane Doe)
Programmable resolution thermometer	23.875 °C
Temperature, Humidity, Barometric Pressure and Lig	391 lux

A red arrow points to the left sidebar containing various icons for document management and analysis.

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Sample name	Location	Status	Pulled on/test date	Pulled by	Reason	Actions
SA_backup_5	Room_1 > Large stability chamber > new Rack > Long_term_plan_2022-L_backup_1	Backup				
SA_backup_6	Room_1 > Large stability chamber > new Rack > Long_term_plan_2022-L_backup_1	Backup				

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Sample name	Location	Status	Pulled on/test date	Pulled by	Reason	Actions
SA_backup_5	Room_1 > Large stability chamber > new Rack > Long_term_plan_2022-L_backup_1	Pulled	24/05/2022 15:58:08	Jane Doe	Need a additional test, the plate breaks.	
SA_backup_6	Room_1 > Large stability chamber > new Rack > Long_term_plan_2022-L_backup_1	Backup				

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ⓘ This plan has been automatically marked as finished on 2021-10-07 12:09:34.
1 test(s) have been marked Executed and 0 test(s) have been marked Not executed.

[Edit](#) [Duplicate](#) [Export PDF](#)

Type: No type Project: reagent Module: Stability Test Status: **Finished**

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Short term stability plan (#6)

Product name:	Product D
Type:	Microbiological
Project:	No project
Module:	Stability test
Condition:	25°C ± 2°C/60% RH ± 5% RH
Status:	Stopped on 25/05/2022 13:43:14 by Jane Doe

Owner:	Jane Doe
Collaborators:	Albert Einstein John Pierce Patricia Perez

Tests				
Day	Stability test	Status	On/by	Sensor(s)
23/05/2022	SA_113	Executed	on 24/05/2022 18:02:54 by Jane Doe	Programmable resolution thermometer: 23.88 °C Temperature, Humidity, Barometric Pressure and Lig: 391 lux
24/05/2022	SA_114	Executed	on 24/05/2022 18:02:59 by Jane Doe	Programmable resolution thermometer: 23.88 °C Temperature, Humidity, Barometric Pressure and Lig: 391 lux
25/05/2022	SA_115	Pending		

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